

Liquid for inhibiting algae growth

Supplied in a 1 liter dosing bottle, **Clinet[®] anti-algae** is used, once diluted to 0.1%, to prevent algae growth in laboratory water baths.

Application

Algae-inhibiting liquid for laboratory water baths.


How to use

Dilute **Clinet[®] anti-algae** to 0.1%. Pour a 20 ml dose into a water bath of 20 liters of distilled water.

Renew the dose every week. Change the water every month.

Presentation

1 liter bottle with dosing dispenser.

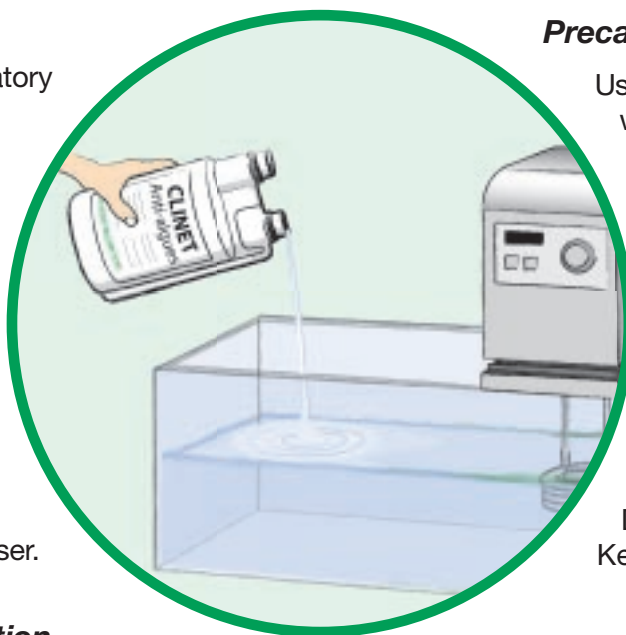

Precautions

Use only demineralized or deionized water, as the algicide activity is reduced by hard water.

Never mix with anion-active detergents and/or strong oxydants (such as bleach...).

In case of contact with the eyes, rinse abundantly with water and consult a doctor.

Do not swallow.
Keep out of reach of children.

Composition

Quaternary ammonium, forms a complex with calcium, preservative.

✂

- Please send me :
- a price offer for Clinet[®] anti-algae.
 - technical information on Clinet[®] anti-algae.
 - a free sample of Clinet[®] anti-algae.

Company

Name

Position

Address

Zip code/City

Country

Tel. Fax

E-mail

Hygiene & Security

information samples

- | | | | |
|-------------------------------|--|--------------------------|--------------------------|
| Clinet[®] 101 | (chlorhexidine cleaning wipes) | <input type="checkbox"/> | <input type="checkbox"/> |
| Clinet[®] 202 | (indivi. chlorhexidine cleaning wipes) | <input type="checkbox"/> | <input type="checkbox"/> |
| Clinet[®] 303 | (bactericidal spray) | <input type="checkbox"/> | <input type="checkbox"/> |
| Anabac[®] | (autoclave deodorant) | <input type="checkbox"/> | <input type="checkbox"/> |
| BioParfum[®] | (odor neutralizers) | <input type="checkbox"/> | <input type="checkbox"/> |

BagSystem[®] for sample preparation

- | | | | |
|------------------------------|------------------------|--------------------------|--------------------------|
| BagMixer[®] | (sample blenders) | <input type="checkbox"/> | |
| BagFilter[®] | (sterile blender bags) | <input type="checkbox"/> | <input type="checkbox"/> |
| BagPipet[®] | (automatic pipette) | <input type="checkbox"/> | |

Distributor

Please fax this request to your local dealer or to :