T.B.X. AGAR

Selective and chromogenic medium for *E. coli* detection in food according to ISO 16649-2.

TYPICAL FORMULA (g/L)

Tryptone	20.0
Bile salts no.3	1.5
Agar	14.0
X-GLUC	75.0 mg
final pH	7.2 +/- 0.2

DESCRIPTION

T.B.X. AGAR is recommended by ISO 16649-2 to detect *Escherichia coli* in products for human or animal consumption by 44 °C incubation and by inclusion or superficial inoculation method. Otherwise, the medium can be inoculated by membrane filtration, according to conventional methods.

PRINCIPLE

Bile salts inhibit the growth of Gram positive bacteria. The chromogenic substratum X-GLUC (5-bromo-4-chlor-3 indolil- β -D-glucuronide), has a differentiating action: the hydrolysis of such substratum, through the β -glucuronidase enzyme, causes a blue-green color pigment formation.

PREPARATION

Suspend 35.6 g powder into 1 liter of distilled or deionized water. Autoclave at 121 °C for 15 min. Cool till about 50 °C. Dispense into Petri dishes.

TECHNIQUE

The procedure recommended by ISO 16649-2 is the following:

- Inoculate into 90 mm plates, 1 ml of sample suspension and 1 ml of its decimal dilutions
- Add 15 ml roughly of medium thermostated at 47-50 °C
- Mix carefully the inoculum with the medium.
- Incubate at 44 °C for 18-24 hours. In case of suspect presence of stressed colonies, pre-incubate for 4 hours at 37 °C before incubating at 44 °C. Do not exceed the temperature of 45 °C.
- Count the blue and green-blue colonies as presumptive *E. coli*
- Indicate results as CFU/g considering the "dilution factor"

INTERPRETATION OF RESULTS

E. coli, among the enterobacteria, is one of the few β -glucuronidase positive species, together with some strains of *Salmonella* and *Shigella*, therefore grows on medium with blue or green-blue colonies. The β -glucuronidase negative enterobacteria grow with colourless colonies.

STORAGE CONDITIONS

10-30 °C away from light, until the expiry date on the label or until signs of deterioration or contamination are evident.

WARNING AND PRECAUTIONS

The product is classified as hazardous by current legislation. It is recommended that the Safety Data Sheets be consulted before use.

The product must be used only by properly trained operators.

Previously to the use be sure the product has not signs of deterioration or contamination.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to current national and local regulations

REFERENCES

ISO 16649-2 (2001): *Microbiology of food and animal feeding stuffs-Horizontal method for the enumeration of E. coli – Part 2: colony count technique at 44°C using 5-bromo-4-chloro-3-indolyl- β -D-glucuronic acid.*

PRODUCT SPECIFICATIONS

NAME

T.B.X. AGAR

PRESENTATION

Dehydrated culture medium.

STORAGE

10-30°C

PACKAGING

Code	Content	Packaging
610224	500 g	500 g powder in plastic bottle
620224	100 g	100 g powder in plastic bottle

pH OF MEDIUM

7.2 ± 0.2

USE

T.B.X. AGAR is recommended by ISO 16649-2 for detecting presumptive *Escherichia coli* in products for human consumption or for animal feeding with incubation at 44 °C and by inclusion or superficial inoculation method.

The medium can be inoculated also by filtration membrane, referring to the conventional methods.

TECHNIQUE

Refer to the technical sheet of the product.

APPEARANCE OF THE MEDIUM

dehydrated medium

Appearance: homogeneous.

Color: creamy white

Ready medium

Appearance: clear

Color: white

VALIDITY

2 years.

QUALITY CONTROL FOR THE USER

1. Control of general characteristics, label and print

2. Sterility control

7 days a 25 ± 1 °C, in aerobiosis

7 days a 36 ± 1 °C, in aerobiosis

3. Microbiological control

Inoculum for productivity: 10-100 UFC/ml

Inoculum for selectivity: 10⁴-10⁵ UFC/ml

Inoculum for specificity: ≤ 10⁴ UFC/ml

Incubation conditions: 44 ± 1 °C for 18-24 hours

Microorganism		Growth	Color of colonies
<i>Escherichia coli</i>	ATCC 25922	good/very good	green-blue
<i>Salmonella typhimurium</i>	ATCC 14028	good/very good	colorless
<i>Enterococcus faecalis</i>	ATCC 19433	Inhibited	

TABLE OF SYMBOLS

SYMBOL	MEANING
	Catalogue number
	Manufacturer
	Kit content
	Use by
	Batch code
	Keep away from heat
	Consult accompanying documents
	Temperature limitation

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